

www.targetingsystems.net

# In Vivo-Targefect

## **Transfection Protocol**

CATALOG NO.	SIZE
IN-VIVO-TARG-001	200 ul
IN-VIVO-TARG-002	600 ul
IN-VIVO-TARG-003	3.2 ml

**Description:** In Vivo-Targefect transfection reagent is a proprietary, biodegradable polymer-based formulation optimized for intramuscular *in vivo* delivery of plasmid DNA and RNA. *The formulation is free from component sof animal origin* 

**Storage:** Store reagent at 4°C upon receipt. If stored properly, reagent is stable for 1 year from date of storage.

## **Recommended Modes of Administration:**

Intramuscular injection

#### **Intended Use:**

For research use only. Not intended for any diagnostic or therapeutic use.

**TABLE 1: DOSING** 

Animal body weight (g)	Transfection Reagent (μl) / per Injection
10 – 35 g (mouse)	20-50 μΙ
35 – 200 g (small rodents)	50-100 μΙ
200 – 450 g (rat)	125-250 µl

### **Recommended Protocol**

Gently mix the In Vivo-Targefect transfection reagent by inverting the tube several times. Use 1 ml conical bottom cryotubes (or Eppendorf tubes) for complex formation.

We recommend trying the following conditions.

**Condition 1:** Add 20 ug of plasmid DNA to each tube, add water to make the volume to 13 ul, mix well by flicking the tube 10-12 times to create a vortex action (do not vortex). Add 7 ul of In vivo-Targefect, mix well again by flicking the tube. Incubate at room temperature for 30 min. Inject 20 ul/mouse intramuscular injection.

**Condition 2:** Add 50 ug of plasmid DNA to each tube, add water to make the volume to 32.5 ul, mix well by flicking the tube 10-12 times to create a vortex action (do not vortex). Add 17.5 ul of In vivo-Targefect, mix well again by flicking the tube. Incubate at room temperature for 30 min. Inject 50 ul/mouse intramuscular injection.

## If working with RNA or siRNA use 50 ug of siRNA in place of DNA and complex as above.

Use Table 1 to determine amount to be injected for larger animals and prepare transfection complexes accordingly.

Assay for target gene expression at desired time intervals. We see target gene expression as early as 2 days following intramuscular. Gene expression typically increases at day 14 and is sustained for at least 1 month in mouse models.

#### **Limited Use Label License:**

The purchase of this product conveys to the purchaser the limited right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. This product is for research purposes only and is not for use in commercial applications of any kind. For information on obtaining additional rights, please contact Targeting Systems at <a href="mailto:info@targetingsystems.net">info@targetingsystems.net</a>

#### **Limited Product Warranty:**

This warranty limits our liability to replacement of this product. No other warranties of any kind, express or implied, including, without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Targeting Systems. Targeting Systems shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product. This product is developed, manufactured, and sold for research purposes only. Reagent is not suitable for administration to humans.